Labortischserie / Lab Bench Series

HPLC

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Contents

1	Overview	15
2	Safety	15
2.1	Important safety instructions	15
2.2	Intended use.....	16
2.3	Non-intended use.....	16
2.4	Examples of foreseeable misuse	16
2.5	Who may use and operate the lifting tables	16
2.6	Safety signs and symbols	17
	Safety symbols on the device and in the user manual.....	17
	Symbols on the nameplate	17
3	Product description	18
3.1	Scope of delivery	18
3.2	Available accessories.....	18
3.3	Model numbers and order codes	18
	Structure of the model number	18
	Structure of the order code	19
	Examples	19
3.4	Technical data	20
	General technical data	20
	Technical data of the height-adjustable version.....	20
4	Handling	21
4.1	Unpacking.....	21
4.2	Transport	21
	Within buildings	21
	Transport over longer distances.....	21
4.3	Commissioning	22
	Positioning	22
4.4	Height adjustment.....	22
	Changing the height.....	22
	Saving and recalling a position	22
	Calibrating the height display / setting an offset	23
	Changing the unit of measurement (centimeter or inch)	23
4.5	Maintenance and care.....	23
	Maintenance	23
	Cleaning	23
4.6	Storage	23
4.7	Final disposal.....	23

1 Overview

- 1 Chemical resistant table top made of TRESPA® TOPLAB® PLUS.
- 2 Control panel for height adjustment and display of table height (height-adjustable models only).
- 3 Feedthroughs for cables and hoses.
- 4 Lifting columns (height-adjustable models) or table columns (models with fixed height).
- 5 Drawer with soft-stop pull-outs (optional).
- 6 Lockable casters.
- 7 Cable duct with integrated socket strip, velcro cable ties and appliance inlet for connecting the mains cable.

2 Safety

2.1 Important safety instructions

- Please read these operating instructions carefully before using the product. It contains important notes on commissioning and handling.
- The user manual is part of the product. Keep it in a safe place. If the product is sold, rented or passed on in any other way, the user manual must be included.
- Heed all warnings. These point out dangers and give recommendations for accident prevention.
- Only install the unit in locations expressly recommended by the manufacturer.
- If condensation forms after unpacking, allow at least 2 hours for the unit to acclimatize.
- This device may only be operated at sockets with protective earth.
- Make sure that you cannot trip over the mains cable and that the cable is not disconnected, is pinched, torn off or squeezed.
- Never reach into the working area of the height adjustment during operation. In addition, make sure that no third person is in the hazardous area during operation. This applies in particular to places that are marked with a corresponding warning symbol. Non-observance can lead to severe crushing or even shearing.
- Only use accessories expressly recommended by the manufacturer.
- Leave all maintenance work to an experienced service technician. Servicing or repairs are required if there is visible damage, if the power cord or plug is damaged, if liquid has been spilled inside the apparatus, if the apparatus has been exposed to rain, if it does not perform as expected, or if it has fallen.
- Only use connection cables or extension cables that are approved for the maximum permissible total current or power.

- Operation of the device in a manner not intended by the manufacturer may affect the protection provided by the device.

2.2 Intended use

- The tables of the HPLC series are products according to the Machinery Directive 2006/42/EC (complete machine).
- The electrically height-adjustable tables of the HPLC series are lifting devices for raising and lowering the integrated work surface.
- The tables are technical work equipment and not intended for private use.
- The table must be placed on a flat, horizontal surface.
- The device may only be used in clean and dry rooms.
- The limit values for temperature and humidity during transport, storage and operation specified under 3.4 „Technical data“ on page 20 must not be exceeded or undercut.
- The height-adjustable table is not suitable for lifting people.
- The table must not be used in potentially explosive atmospheres, especially in areas enriched with oxygen, combustible anesthetics (class AP) or combustible anesthetics with oxidizing agents (class APG).
- Use the table only within the specifications specified in this manual.
- The valid accident guarding regulations according to each country must be considered.
- The device may only be repaired by a specialist.

2.3 Non-intended use

- A non-intended use is given when the equipment is treated in any way that is contrary to any of the points from the chapter 2.2 „Intended use“ on page 16.
- If the lifting tables are not used as intended, handled improperly and used, assembled or handled by untrained personnel, hazards may arise for personnel from the lifting tables.
- The moving of persons with this lifting table, as an example of improper use, is prohibited. In case of improper use the liability of the manufacturer as well as the general operating permit of the lifting table expires.

2.4 Examples of foreseeable misuse

- Overloading of the table by exceeding the maximum permissible mass.
- Outdoor use.
- Use in environments with high humidity > dew point.
- Use in rooms with explosive atmospheres according to ATEX directive.
- Opening the lifting columns or the electrical system of the table.
- Use on insufficiently safe ground.
- Use with damaged supply lines or mechanics.

2.5 Who may use and operate the lifting tables

- Physically and mentally healthy persons who have fully read and understood the instructions for use may use and operate these tables as long as they are not affected by medication or drugs.

2.6 Safety signs and symbols

Safety symbols on the device and in the user manual

The “general mandatory sign” indicates that you should behave attentively. Marked information in this user manual is intended to draw your particular attention. You will receive important information on functions, settings and procedures. Failure to do so may result in malfunction of the equipment or its surroundings.

The “general warning sign” indicates danger points or possible dangerous situations. Marked information in this user manual applies to your special attention. You will receive important information on safety-relevant matters. Failure to do so may result in personal injury, malfunction of the device or the environment.

The warning sign “Hand Injury Warning” indicates that hands or other parts of the body may be crushed, pulled in or otherwise injured at the points marked. It is essential that you observe the safety instructions in the operating instructions.

The warning sign “Warning of electrical voltage” indicates potential hazards due to electrical voltage. The positions on the device marked with this symbol and the information in these operating instructions are intended to draw your particular attention. You will receive important information on safety-relevant matters. Failure to do so may result in personal injury, malfunction of the device or the surrounding area.

Symbols on the nameplate

Conformity to low voltage and EMC directives

Do not dispose of in household waste

Indoor use only

Appliance class I

3 Product description

3.1 Scope of delivery

- 1 x Laboratory table of the HPLC series. Model and accessories according to model number and order code.
- 1 x Region-dependent power cable (see 3.4 „Technical data“ on page 20).
- 1 x This user manual.

3.2 Available accessories

Drawer module “D1”

Standard drawer module. Drawer height: 68mm, front panel height: 119mm.
The width and depth of the drawer depends on the table dimensions.
Loading capacity: 15kg

Waste bottle tray “WT”

Tray with removable anti-slip mat, e.g. for waste bottles which moves along with the tabletop. Depth: 23cm Width: Table width - 20cm, tub height: 5cm.
The width of the tray depends on the table dimensions.

Fixed floor plate “FP”

Chemical resistant shelf made of TRESPA® TOPLAB® PLUS. The dimensions of the floor plate depend on the table dimensions.

3.3 Model numbers and order codes

Structure of the model number

The model number consists of the essential features of the table. Properties in brackets are optional. The structure is as follows:

Series **Width** **Depth** . **Height** (**Travel**) **Region** (. **Power strip**)

Series	“H” as identification of the HPLC product series.
Width	Three-digit width of the table in centimeters. Available widths: 60cm (stocked) 80cm (on request) 100cm (on request) 120cm (stocked)
Depth	Two- or three-digit depth of the table in centimeters. Available Depths: 70cm (on request) 80cm (stocked) 90cm (on request) 100cm (on request)

Height / (travel)	Two-digit basic height and travel of the table in centimeters. For the height-adjustable models, the lowest possible height is indicated as "Height". For models with a fixed height, "Travel" is not specified.												
	<table> <thead> <tr> <th>Available heights (fixed-height models):</th> <th>Available heights (height-adjustable models):</th> </tr> </thead> <tbody> <tr> <td>60cm (stocked)</td> <td>53-83cm (stocked)</td> </tr> <tr> <td>70cm (on request)</td> <td>58-93cm (on request)</td> </tr> <tr> <td>80cm (stocked)</td> <td>63-103cm (stocked)</td> </tr> <tr> <td>90cm (on request)</td> <td>68-113cm (on request)</td> </tr> <tr> <td>100cm (on request)</td> <td>73-123cm (on request)</td> </tr> </tbody> </table>	Available heights (fixed-height models):	Available heights (height-adjustable models):	60cm (stocked)	53-83cm (stocked)	70cm (on request)	58-93cm (on request)	80cm (stocked)	63-103cm (stocked)	90cm (on request)	68-113cm (on request)	100cm (on request)	73-123cm (on request)
Available heights (fixed-height models):	Available heights (height-adjustable models):												
60cm (stocked)	53-83cm (stocked)												
70cm (on request)	58-93cm (on request)												
80cm (stocked)	63-103cm (stocked)												
90cm (on request)	68-113cm (on request)												
100cm (on request)	73-123cm (on request)												
Region	Two-digit regional code. Possible values: "EU", "CH", "UK" and "US". Determines the type of power cable supplied and the type of power supply. For more information see 3.4 „Technical data“ on page 20.												
(Power strip)	Type of the integrated power strip. If the specification is missing, the value specified under "Standard" applies. Standard: 12 x IEC-60320 C13 S1: 9 x Schuko (CEE 7/3) (DE) S2: 8 x T23 (SEV 1011) (CH) S3: 7 x BS1363 (UK) S4: 12 x NEMA 5-15 (USA)												

Structure of the order code

The order code consists of the model number of the table and - separated by hyphens - the given codes of the included accessories.

Model number - **Accessory 1** - **Accessory 2** - ... - **Accessory n**

Examples

Model number: H06080.5330EU | Order code: H06080.5330EU-D1

HPLC laboratory table, width: 60cm, depth: 80cm, base height: 53cm, travel: 30cm, region: EU, drawer module "D1"

Model number: H12080.80CH.S3 | Order code: H12080.80CH.S3-D1-FP

HPLC laboratory table, width: 120cm, depth: 80cm, fixed height: 80cm, region: CH, power strip S3, drawer module "D1" + fixed floor plate "FP"

3.4 Technical data

General technical data

Dimensions	See model number	
Weight	80-140kg depending on version	
Load capacity	250kg with even load distribution 150kg with one-sided load distribution	
Permissible ambient conditions	Temperature	Humidity
Transport	-20 - 60°C / -4 - 140°F	5-90% (non-condensing)
Storage	10 - 40°C / 50 - 104°F	10-80% (non-condensing)
Operation	10 - 30°C / 50 - 86°F	10-80% (non-condensing)
Power cable		
EU	CEE 7/4 plug (Schuko plug) to IEC 60320 C13, length: 5m	
CH	T12 plug (according to SEV 1011:2009) to IEC 60320 C13, length: 5m	
UK	Type G plug (BS 1363) to IEC 60320 C13, length: 5m	
US	NEMA-5 plug to IEC 60320 C13, length: 5m	
Supply voltage		
EU, CH, UK	207-253V / 50Hz	
US	90-127V / 50-60Hz	
Power rating		
Combined power input	2500W max.	
Appliance class	Class I (earth conductor)	

Technical data of the height-adjustable version


Power rating	
Standby	<0,6W
Height adjustment	288W max.
Maximum lifting speed	9mm/s
Duty cycle	10% (S3) -> 1 minute operation per 10 minutes

4 Handling

4.1 Unpacking

Make sure that there are at least two of you when unpacking the table to ensure that the table cannot tip over or fall down the side of the ramp. Depending on the model, the table may exceed a weight that can be safely held by one person alone.

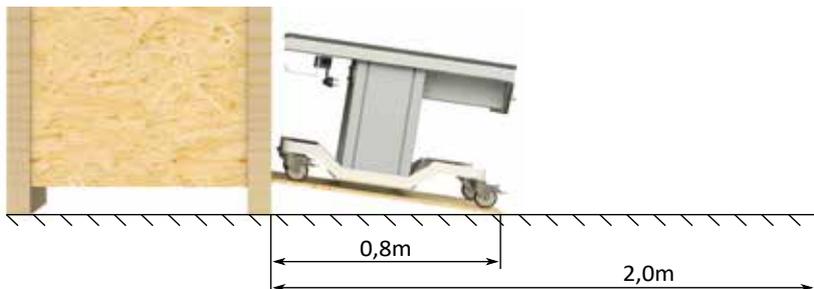If the table is delivered in a wooden box, please observe the following instructions.

The wooden box is equipped with an integrated loading ramp. Position the transport box on an even surface so that there is sufficient space to fold down the ramp and roll out the table.

Open the box lock and fold down the loading wall. Release the parking brakes on the casters and carefully move the table down the ramp.

4.2 Transport

Within buildings

Make sure that all wheel brakes are released and that the connecting cable is unplugged or long enough.

Especially with higher set-ups and / or with extended lifting columns, the table can become unstable or even tip over during transport. If possible, move the table only in the lowest possible position and, if necessary, secure the table and / or the equipment on top of the table against tilting.

The table has a low ground clearance of approx. 2cm in order to be able to occupy the lowest possible position. When moving over for example a door thresholds, it may be necessary to lift the table slightly to avoid sitting on it.

Transport over longer distances

- When transporting over longer distances, especially in vehicles, make sure that the table is secured against transport damage. It is best to use the original packaging.
- The table may only be transported over longer distances in the lowest possible position.
- If the table is equipped with moving parts, such as drawers, these must be secured against slipping.

4.3 Commissioning

Positioning

Position the table in the designated place. With the height-adjustable models, make sure that there is enough space upwards even when the table is fully extended. If possible, fix the casters to prevent unintentional movement.

When connecting the power cable, make sure that it is of sufficient length. If necessary, use an extension cable. Lay the cable in such a way that it cannot be bent or crushed. Also make sure that the cable is laid in such a way that no one can trip over it.

Connect the table to a suitable wall outlet using the supplied power cable. The appliance inlet is located in the cable tray above the integrated power strip.

4.4 Height adjustment


Changing the height

Never reach into the working area of the height adjustment during operation. In addition, make sure that no third person is in the danger area during operation. This applies in particular to places which are provided with a corresponding warning symbol. Non-observance can lead to severe crushing or even shearing.

Make sure that there are no objects in the adjustment range. The table or any objects colliding with it may be damaged or destroyed. Also make sure that the power cable to the table is long enough to fully adjust the table.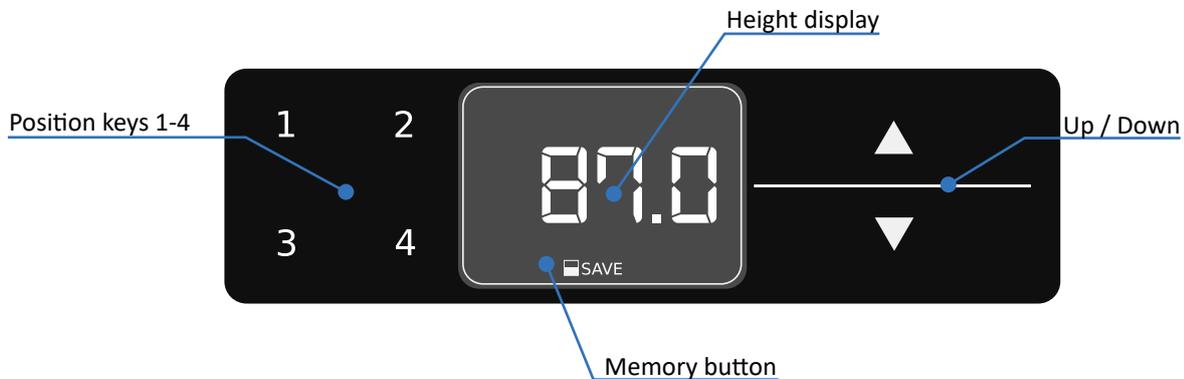

With the electrically height-adjustable versions, the table height can be adjusted via the control panel located under the table top. The display in the middle shows the current table height in cm or inch. Use the two arrow keys to the right of the display to move the table up or down.

If "HOT" is shown in the display, the motor is overloaded due to too long travel. In this case, wait a few minutes until the message disappears automatically.

Saving and recalling a position

The position keys 1 - 4 can be used to move to programmed table heights or to save the current table height. To save the currently set position, press the save key. The display will now show "S-". Press the desired position key (1-4). The controller will now click twice and the display will show, for example, "S1". To move to one of the stored positions, hold down the corresponding position key 1, 2, 3 or 4 until the stored position is reached.

Calibrating the height display / setting an offset

If the display should show a relative height instead of the absolute height, an offset can be set. To change the height shown on the display, press the Memory button. The display will show "S -". Press the Down button for about 5 seconds. The display flashes. You can now use the "Up" and "Down" key to set the desired value. Press the Memory button again to save the currently displayed relative height.

Changing the unit of measurement (centimeter or inch)

Press the position keys 1 and 2 and the "Up" key at the same time. The display will now show "S" and a number (e.g. "S 7"). Press the Up button until "S 5" appears on the display. Press the Save button to confirm the setting. If the display was previously set to centimeter, it will now change to inches or vice versa.

4.5 Maintenance and care

Maintenance

The product is basically maintenance-free but not wear-free. If a wear-related defect should occur, please contact your dealer or Sonation directly to clarify the further steps.

Cleaning

Use a soft, damp cloth to clean the surfaces. Aggressive solvents, bases or acids must not be used for cleaning.

4.6 Storage

- Height-adjustable models should be moved to the lowest possible position for storage.
- Disconnect the table from the mains.
- Secure moving parts such as drawers to prevent accidental extension.
- If possible, use the original packaging for storage.
- The ambient conditions listed under 3.4 „Technical data“ on page 20 must be observed.

4.7 Final disposal

Dispose the table in accordance with local and national safety and environmental regulations.

The table contains electronic components, cables, metals, plastics, etc. and has to be disposed in accordance with applicable environmental regulations of the respective country. The disposal of the product in Germany is subject to the Elektro-G (RoHS) and in the European area of the EU Directive 2002/95/EC or the respective national legislation.

The table can also be returned to the seller for disposal. In this case the cost of transport, including the packaging costs shall be borne by the buyer. Please notify the seller of the return within 14 days before the expected arrival.

IMPROVE YOUR SCIENCE

Sonation GmbH
Alte Schulstr. 39
88400 Biberach
Germany