USBB-S
USBB-M
USBB-L
USBB-XL

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according to 2006/42/EC and 2004/108/EC

Manufacturer:
Sonation GmbH
Alte Schulstr. 39
88400 Biberach
Germany

Mr. Christof Voelkle is authorized to compile the technical file.
Address: Sonation GmbH, Alte Schulstr. 39, 88400 Biberach

Hereby we declare that the following noise reduction boxes

USBB-S, USBB-M, USBB-L, USBB-XL

comply with all relevant provisions of low voltage directive
2006/95/EC and EMC directive 2004/108/EC.

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CONTENT

Important safety instructions.....	8
Operating conditions and conventional usage.....	8
Models.....	9
Technical data.....	9
Scope of supply.....	9
Accessories.....	9
Overview.....	10
Installation.....	10
Operation.....	10
Maintenance.....	10
Notes.....	11

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NOTIZEN

A large grid of graph paper for taking notes, consisting of 20 columns and 30 rows of small squares. The word 'NOTIZEN' is printed in a bold, black, sans-serif font in the top-left corner of the grid, within the first row and the first four columns.

IMPORTANT SAFETY INSTRUCTIONS

- Read these instructions.
- Keep these instructions.
- Heed all warnings.
- Do not block any ventilation openings. Install in accordance with the manufacturer's instructions.
- Use only with accessories specified by the manufacturer, or sold with the apparatus.
- Refer all servicing to qualified service personnel. Servicing is required when the apparatus has been damaged in any way, such as power-supply cord or plug is damaged, liquid has been spilled or objects have fallen into the apparatus, the apparatus has been exposed to rain or moisture, does not operate normally, or has been dropped.
- If the apparatus is used in a manner not specified by the manufacturer the protection provided by the equipment may be impaired.

OPERATING CONDITIONS AND CONVENTIONAL USAGE

- The box is exclusively designed as a noise reduction device for ultrasonic baths.
- The ultrasonic bath may not have tangible dangerous parts.
- The box must be installed on an even and horizontal surface.
- The exhausted airflow must be able to stream unhindered. The rear air discharge opening must be distant from other objects at least 15 cm.
- The fan must be able to turn free and may not be blocked.
- The box may only be used in clean and dry areas.
- During transport and storage temperature and humidity must be within the following range:
Transport: Temperature: -20 to 60°C (-4 to 140°F), humidity: 5 to 90%
Storage: Temperature: 10 to 40°C (50 to 104°F), humidity: 10 to 80%
- The box may not be operated above 30°C / 86°F ambient temperature and above 80% humidity.
- The inner temperature of the box may not exceed 50°C / 122°F
- If condensation water is formed after unpacking the box, an acclimatisation period of 2 hours must be adhered prior to installation.
- The box must be kept away from easily inflammable and combustible liquids.
- If a liquid should penetrate the electrical part of the box it is to be set immediately out of operation and must be checked by a specialist.
- The valid accident guarding regulations according to each country must be considered.
- A repair of the box may be accomplished only by a specialist.

MODELS

Model	Inner dims WxDxH	Outer dims WxDxH	Weight
USBB-S:	454x269x316 mm	512x350x425 mm	15kg
USBB-M:	479x329x415 mm	557x410x545 mm	20kg
USBB-L:	529x379x496 mm	607x460x633 mm	25kg
USBB-XL:	782x610x671 mm	704x529x516 mm	30kg

TECHNICAL DATA

Voltag / Current Consumption:

12V DC, 1000mA max

Disconnecting device:

The wall power supply serves as the disconnecting device. To completely disconnect the equipment from the electric supply network disconnect the wall power supply. Please assure that the wall power supply is accessible at all times.

Sound proofing:

Environmental friendly recycling product. Oil and water repellent. Inflammability fulfills FMVSS 302 and DIN 75200.

SCOPE OF SUPPLY

- One noise reduction box USBB-S, USBB-M, USBB-L or USBB-XL
- One wall power supply 12V, 1000mA

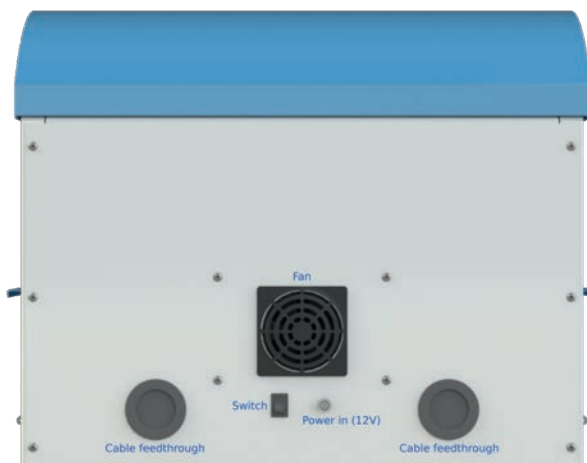
ACCESSORIES (NOT INCLUDED)

Side table for easy accessibility

Especially with the larger models USBB-L and USBB-XL accessibility for smaller people is difficult when the system is set up on a table with standard height. For this purpose we offer a lower side table with the dimensions 610x800x620mm (WxDxH).

Order code: USBB-Easy table

OVERVIEW

INSTALLATION

- Place the noise reduction box at the position where the ultrasonic bath will be operated. Plug the wall power supply into a wall outlet and the plug on the cord to the jack on the back of the noise reduction box.
- Open the noise reduction box by pushing the upper handle backwards and position the ultrasonic bath inside the noise reduction box.
- Feed the power cord of the ultrasonic bath through one of the two rubber feedthroughs on the back
- Make sure that the fan is working. To do so turn on the switch on the back panel. Please make sure that the fan is not blocked and is at least 5cm (2") away from the wall or other equipment. The noise reduction box is now ready to use and the ultrasonic bath can be powered and operated.

OPERATION

Use your ultrasonic bath as described in the users manual. If the heater of the ultrasonic bath is used, the fan must be turned on.

MAINTENANCE

To clean the surface please use a soft and moist tissue.

The function of the fan should be checked every month. Also the fan guard must be checked if dust or other sorts of pollution could block the air flow. Depending on ambient conditions, this may be necessary earlier or.

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IMPORTANT NOTICE TO CONSUMERS

This product has been manufactured according to strict specifications and voltage requirements that are applicable in the country in which it is intended that this product should be used. If you have purchased this product via the internet, through mail order, and/or via a telephone sale, you must verify that this product is intended to be used in the country in which you reside.

WARNING: Use of this product in any country other than that for which it is intended could be dangerous and could invalidate the manufacturer's or distributor's warranty. Please also retain your receipt as proof of purchase otherwise your product may be disqualified from the manufacturer's or distributor's warranty.